

"Because Pedigree Matters"



Laboratory **Filter Media**



In technical collaboration with
Hahnemühle 
Life Science

Borosil is India's
most trusted laboratory brand



SINCE

1962

Borosil is born as a Public Limited Company. The company begins to sell products under the "Borosil" brand.

1963

Corning Glass Works, USA joins Borosil Ltd as a Technical and Financial Collaborator.

2008

Vyline Glassworks sets up a state of the art manufacturing facility for laboratory and consumer glassware. in Bharuch, Gujarat

2013

Vyline sets up a second manufacturing unit for laboratory glassware in Tarapur, Maharashtra.

2015

Borosil launches Labquest - its entry into laboratory instruments.

2016

Borosil acquires controlling interest in Klasspack Pvt Ltd, in Nasik, Maharashtra and enters the pharmaceutical primary packaging market.

2017

Gujarat Borosil produces the world's first fully 2 mm solar glass, facilitating a revolutionary 30% increase in solar module efficiency.

2019

Borosil Technologies Limited with a mandate to research, design, develop laboratory equipment is set up at Pune, India under the brand Labquest.



We are very excited to announce our collaboration with Hahnemühle Fine Art Germany, pioneers of filter papers since 1883.

There are very few companies worldwide that can be proud of such a long history like Hahnemühle.

Borosil's vision is to be the most customer centric company in India.

In keeping with this vision, this strategic partnership is to provide Indian scientists and upcoming intellects with quality filter papers that will aid in producing accurate and consistent results.

Shreevar Kheruka
Managing Director

BOROSIL LIMITED

- 1584** Establishment of Hahnemühle
- 1883** Filter papers produced for the first time
- 1927** TO **2004** Hahnemühle was part of the 'Schleicher & Schuell' group
- 2008** Pure filter papers are marketed directly under the Hahnemühle name



One of the pioneers in the filter paper industry, Hahnemühle, a company headquartered in Dassel, Germany has been in the filter paper business for more than 130 years. The papers, since then, have been used in a wide range of industries including the food and beverage industry, the pharmaceutical and chemical industry, and agriculture, environmental monitoring and automotive engineering.

The strict QC, stringent parameters and paper specifications that have been persisting since decades, have made the papers gain a reputation of producing reproducible filtration results, thanks to the high batch to batch uniformity.

Hahnemühle has been an official "Brand of the Century" since 2016, and therefore belongs to the exclusive circle of Germany's strongest brands.



We are very excited about our new strategic partnership with Borosil for the APJ market, specifically India, initially.

With Borosil, we have met a very committed team here that is ahead-of-the-curve in the Life Science market, committed, competent and single-minded.

Hence, we together pursue this growth opportunity for both of us, establish high-quality filter papers "Made in Germany" in the Indian market.

We have great confidence in the entire operational and leadership team that together we will provide significant value to the Borosil customers in India to further differentiate Borosil in the Indian market.

Jan Wölfle
CEO

Hahnemühle FineArt GmbH

The benchmark in Filter Papers



Borosil filter papers are made in collaboration with **Hahnemühle** - the pioneers in speciality papers since **1883!**

The papers contain upto **98% alpha cellulose**, the most stable form of cellulose in paper pulp, hence making the papers highly **stable and durable**. To ensure batch to batch **uniformity and consistency** in filtering performance, no recycled materials are used in the manufacturing processes. Very **low ash content**(Lowest value 0.004%) aid the laboratory professionals in accurate quantitative analysis with minimal errors. **Made in Germany**, the manufacturing processes and product specifications are compliant to the stringent **European QC** standards, guaranteeing the users finest quality filter papers.



Content

High purity Qualitative filter papers

For qualitative analysis in laboratories

6 - 9

Ashless Quantitative filter papers

Suited for routine routine gravimetric tests and sample preparation for instrumental analyses.

10 - 13

Hardened Quantitative filter papers

Specially recommended for filtration applications requiring vacuum and pressure filtration, and for the use of acidic and alkaline solutions.

14 - 17

Glass Fibre Filters

100% micro-borosilicate glass fibres for specific application in environmental analyses

18 - 19

Thimbles

Cellulose and Glass Fiber thimbles for accurate and consistent extraction applications.

20 - 22

Industrial Applications

24 - 37



Qualitative

Filter Papers

Highly pure filter paper
(approx. 0.08% ash)
recommended for
qualitative analyses in laboratories.

Made of super- refined cotton linters and cellulose.

Alpha-Cellulose content upto 98%, therefore ensuring high stability and durability of the papers

Very low content of minerals (approx. 0.08% ash)

Ideal for precise identification of materials and for sample preparation prior to sensitive detection methods

Specifications

Grade	Sample	Speed	Retention range[μm]*	Filtration Herzberg (s)	Thickness (mm)	Weight [g/m ²]
B59501	Medium fine precipitates	Medium fast, thin	4 – 7	160	0.15	68
B59702	Medium-fine, crystalline precipitates	Medium fast	4 – 7	155	0.18	85
B59303	Fine precipitates	Medium slow	2 – 5	450	0.17	84
B60404	Coarse precipitates	Fast	12 – 25	50	0.19	79

* Approximate values, ** Measured with 100mm water column instead of 50 mm

Applications

Grade B59501

- For medium-fine, crystalline precipitates
- For determining the overall fat content of food products
- Determination of the unsaponifiable fraction in fats and oils
- Digestion of solids with aqua regia
- Sample preparation for analysis of beer/malt and beer based beverages.
- Particle separation in Juices, Wine
- Qualitative analysis of milk and milk products.
- Sample preparation for water analysis

Grade B59303

- For fine crystalline precipitates
- Identification of Barium sulphate, tin sulphide
- Soil analysis
- Particle separation in Juices, Wine

Grade B59702

- For medium-fine, crystalline precipitates, Calcium oxalate, metal sulphide
- Determination of the coagulable proteins in beer, malt industries
- Determination of the grade of fermentation in beer, malt, beer based beverages.
- Determination of the unsaponifiable fraction in fats in edible oils and fat industries.
- Determination of the fat content of food, milk and dairy products
- Preparation of samples and removal of CO₂ in the beverage industry
- Particle separation in Juices, Wine

Grade B60404

- For coarse crystalline precipitates
- Sodium chloride in foodstuffs, ferrous hydroxide, aluminium hydroxide and metal sulphide analysis
- Routine cleaning of organic extracts and biological fluids
- Qualitative analysis of milk and milk products
- High flow rates in air pollution monitoring and exhaust fumes detection
- Particle separation in Juices, Wine and other beverages



Ordering Information

Grade name	Size (dia in mm)	Pack size	Product code
B59501	70	100	FPB595010070
B59501	90	100	FPB595010090
B59501	110	100	FPB595010110
B59501	125	100	FPB595010125
B59501	150	100	FPB595010150
B59501	185	100	FPB595010185
B59501	240	100	FPB595010240
B59702	70	100	FPB597020070
B59702	90	100	FPB597020090
B59702	110	100	FPB597020110
B59702	125	100	FPB597020125
B59702	150	100	FPB597020150
B59303	70	100	FPB593030070
B59303	90	100	FPB593030090
B59303	110	100	FPB593030110
B59303	125	100	FPB593030125
B59303	150	100	FPB593030150
B60404	47	100	FPB604040047
B60404	55	100	FPB604040055
B60404	70	100	FPB604040070
B60404	90	100	FPB604040090
B60404	110	100	FPB604040110
B60404	125	100	FPB604040125
B60404	150	100	FPB604040150
B60404	185	100	FPB604040185
B60404	240	100	FPB604040240
B60404	320	100	FPB604040320

Filter paper SHEETS

Grade name	Size (in cm)	Pack size	Product code
B59501	46*57	100	FPB595014657
B59702	46*57	100	FPB597024657
B59303	46*57	100	FPB593034657
B60404	46*57	100	FPB604044657



Ashless Quantitative

Filter Papers

Ashless filters (avg ash content 0.004 %) for quantitative analysis, routine gravimetric tests and sample preparation for instrumental analyses.

- Made of super - refined cotton linters and cellulose
- Average ash content 0.004%
- Acid-washed and rinsed with water to neutralise it
- Free of minerals and metallic ions,hence ideal for the detection of metallic ions

Specifications

Grade	Sample	Speed	Retention range[μm]*	Filtration Herzberg (s)	Thickness (mm)	Weight [g/m ²]
B58940	Fine crystalline precipitates	Medium fast, thin	2 - 4	450	0.17	84
B58941	Coarse and gelatinous precipitates	Medium fast	12 - 25	50	0.19	79
B58942	Very fine crystalline precipitates	Medium slow	<2	750**	0.16	84
B58943	Coarse precipitates	Fast	4 - 12	140	0.18	85
B58944	Fine crystalline precipitates	Slow thin	2	900	0.15	74

* Approximate values, ** Measured with 100mm water column instead of 50 mm

Applications

Grade B58940



- For fine crystalline precipitates
- Determination of sulfates, carbonates and organic materials
- Analysis of oil/fats in edible oil and fat industries.
- Measurement of trace elements in soil and fertilizers.
- Measurement of nitrogen in soil and fertilizers testing.
- Determination of free amino acids and total amino acids in soil and fertilizers.
- Determination of K and P in soil and fertilizers
- Gravimetric analysis of meat and meat products, chemicals, pharmaceuticals, animal feed and in Oil refineries.
- Measurement of non metallic inorganic compounds in water analysis

Grade B58941



- Fast filtration for coarse and gelatinous precipitates.
- Gravimetric analysis in sugar, milk and milk products, meat and meat products, animal feed products, chemicals
- Total dry/ash residue in samples
- For food analysis
- Blaine test in cement
- Analysis of beverages.
- Preparing fruit juice samples for photometric measurements (e.g. phosphate)
- Determination of solids and turbidity (Feld method) in beer, malt, beer based beverages.
- Determination of proteins in wort and beer via $MgSO_4$ precipitation.
- Analysis of ash content in foodstuffs.
- Measurement of trace elements in soil and fertilizers.
- Determination of K and P in soil and fertilizers.
- Measurement of non metallic inorganic compounds in water analysis

Grade B58942



- For very fine crystalline precipitates
- Analyses of oil/fats: proportion of soluble contaminants
- Determination of chemical elements, radioactive trace elements in water used for the beverage industry.
- Measurement of trace elements, free amino acids, total amino acid in soil and fertilizers.
- Measurement of soluble sulphate in soil and fertilizers.
- Determination of K and P in soil and fertilizers
- Measurement of non metallic inorganic compounds in water analysis
- Gravimetric analysis of chemicals, pharmaceuticals, milk and milk products, , eat and meat products

B58943



- Medium-fast filtration for coarse precipitates
- For food analysis
- Determination of proteins in wort and beer via $MgSO_4$ precipitation
- Analysis of alkaline earth carbonates and galvanic baths
- Blaine test of cement
- Analysis of beverages: determination of nitrogen compounds by phosphor molybdenum and carbohydrates by hydrolysis
- Measurement of trace elements, free amino acids, total amino acid in soil and fertilizers.
- Determination of K and P in soil and fertilizers
- Gravimetric analysis of animal feed, chemicals, pharmaceuticals, milk and milk products, sugar, meat and meat products

Grade B58944



- For fine crystalline precipitates.
- Analysis of CaC_2O_4 , $PbSO_4$, $BaSO_4$ (hot-felled precipitates)
- Measurement of trace elements in soil and fertilizers.
- Determination of K and P in soil and fertilizers
- Gravimetric analysis of chemicals, pharmaceuticals, animal feed

Ordering Information

Grade name	Size (dia in mm)	Pack size	Product code
B58940	70	100	FPB589400070
B58940	90	100	FPB589400090
B58940	110	100	FPB589400110
B58940	125	100	FPB589400125
B58940	150	100	FPB589400150
B58940	185	100	FPB589400185
B58940	240	100	FPB589400240
B58941	70	100	FPB589410070
B58941	90	100	FPB589410090
B58941	110	100	FPB589410110
B58941	125	100	FPB589410125
B58941	150	100	FPB589410150
B58941	185	100	FPB589410185
B58942	55	100	FPB589420055
B58942	70	100	FPB589420070
B58942	90	100	FPB589420090
B58942	110	100	FPB589420110
B58942	125	100	FPB589420125
B58943	12.7	1000	FPB589430012
B58943	125	100	FPB589430125
B58944	90	100	FPB589440090
B58944	110	100	FPB589440110
B58944	125	100	FPB589440125
B58944	150	100	FPB589440150
B58944	185	100	FPB589440185

Filter paper SHEETS

Grade name	Size (in cm)	Pack size	Product code
B58940	46*57	100	FPB589404657
B58941	46*57	100	FPB589414657



Hardened Ashless

Filter Papers for quantitative analysis :

Hardened ashless filters (ash content approx. 0.003%), especially recommended for filtration applications requiring vacuum and pressure filtration, and for the use of acidic and alkaline solutions.



In technical collaboration with
Hahnemühle



- Made of super-refined cotton linters and cellulose
- Acid-washed and free of minerals and metallic ions, ideal for the detection of metallic ions
- Very high wet strength; easy to scrap off or wash off precipitates
- Hardened by small quantities of nitrogen-containing resin, which does not introduce significant impurities into the filtrate
- High resistance to aggressive chemical components, like sulphuric and nitric acids (up to 40 % at 50 ° Celsius) and alkalis (up to 10 % at 20 ° Celsius)
- For analytical applications, routine quantitative and / or gravimetric procedures

Specifications

Grade	Sample	Speed	Retention range [μm]*	Filtration Herzberg (s)	Thickness (mm)	Weight [g/m^2]
B15540	Fine crystalline precipitates	Medium fast, thin	4 - 12	170	0.16	90
B15541	Coarse crystalline precipitates	Fast	12 - 25	50	0.17	88
B15542	Very fine crystalline precipitates	Slow	≤ 2	600**	0.14	90

* Approximate values, ** Measured with 100mm water column instead of 50 mm

Applications

Grade B15540

- For fine crystalline precipitates
- Gravimetric determination of metals in acidic/alkaline solutions, quality control of chemicals.
- Gravimetric analysis in oil refinery

Grade B15542

- For very fine crystalline precipitates
- Gravimetric analyses of fine metals: barium and lead sulphate, nickel and tin sulphide, oxalate and calcium fluoride
- Gravimetric analysis in oil refinery

Grade B15541

- For coarse crystalline precipitates
- For gravimetric tests in cement, mining industry
- Analysis in electroplating: baths of aluminium, chrome and copper



Ordering Information

Grade name	Size (dia in mm)	Pack size	Product code
B15540	125	100	FPB155400125
B15540	150	100	FPB155400150
B15541	125	100	FPB155410125
B15542	125	100	FPB155420125
B15542	150	100	FPB155420150





Glass Fiber

Filters

100% micro borosilicate glass fibres for specific applications
in environmental analysis



In technical collaboration with
Hahnemühle



- Made of 100% micro-borosilicate glass fibres
- Chemically stable in acidic solutions (except hydrofluoric acid) and alkaline solutions in moderate concentrations
Extremely low metal content
- Maintains its properties up to 500°C
- High flow speed and high permeability to air

Specifications

Type	Air permeability* (Resistance** mbar) [L/m ² s]	Retention rate % NaCl-particle size <1µm***	Filtration Gurley [s]	Weight [g/m ²]	Thickness (mm)	Max T[°C]	Retention range[µm]#
BGF50A	25	99.97	19	56	0.29	500	1-3
BGF52C	54**	99.995	25	54	0.28	500	1-3

* as per DIN 53887

** Air resistance at 400cm³/s, A = 10cm²

*** Tested with NaCl particles size <1 µm, main fraction at 0.3 to 0.5 µm

#Retention range are approximate values.

BOROSIL[®] Scientific Laboratory Filter Media

Thimbles

Cellulose and Glass Fiber thimbles for accurate and consistent extraction applications.



Cellulose and Glass fibre thimbles for extraction of components

- Extraction thimbles to be used in your regular soxhlet apparatus based applications
- Consistent, high porosity of thimbles ensuring a rapid flow rate
- Made of cellulose or glass fibre based on application

Applications - Cellulose thimbles

- Extraction of fatty, greasy materials in foodstuffs, paints, varnishes etc
- Extraction of active agents from pharmaceuticals and plastic softening agents
- Analysis of pesticide waste, polyaromatic carbohydrates and dioxins in foodstuffs.
- Determination of oil content in oil-bearing seeds

Applications - Glass Fibre thimbles

- Extraction of solvents which are incompatible with cellulose fibres
- Gravimetric collection of dust particles or aerosols from hot air and gas flows
- Extraction during special biochemical analysis

Technical Data

Product	Material	Wall Thickness (mm)	Weight (gsm)	Air Permeability (L/m ² sec)
Cellulose Thimble	100% cellulose without binder	1.6	2.7	25
Glass Fibre Thimble	100% Micro-glass fibre without binder	1.3	3	15

Ordering information

Product Codes	Description	Size (Inner dia x height)	HSN Code	Pack size
FPBCTH01990	BCTH 19x90mm Cellulose Thimbles	19x90	48237090	25
FPBCTH022080	BCTH 22x80mm Cellulose Thimbles	22x80	48237090	25
FPBCTH022100	BCTH 22x100mm Cellulose Thimbles	22x100	48237090	25
FPBCTH025080	BCTH 25x80mm Cellulose Thimbles	25x80	48237090	25
FPBCTH025100	BCTH 25x100mm Cellulose Thimbles	25x100	48237090	25
FPBCTH028100	BCTH 28x100mm Cellulose Thimbles	28x100	48237090	25
FPBCTH033080	BCTH 33x80mm Cellulose Thimbles	33x80	48237090	25
FPBCTH033100	BCTH 33x100mm Cellulose Thimbles	33x100	48237090	25
FPBGFTH01990	BGFTH 19*90mm Glass Fibre Thimbles	19x90	70197100	25

Air Monitoring Filters

The increasing presence of particulate matter and pollutants poses severe risks to human health and the environment.

Protecting Earth's vital resource, air, is a global priority.

- Designed to detect and measure pollutants such as PM2.5, PM10, sulfur dioxide, fluorides, heavy metals, and radioactive substances.
- Utilizes gravimetric analysis for precise measurement.
- Employs membrane filter discs and filter holders for effective monitoring.

Ordering information

Product Codes	Size	Pack size
FPBGFEP2000	203*254mm sheet (8*10inches)	100
FPBPTPM25	PTFE PM2.5 filter discs 2 µm pore size, 46.2 mm dia.	100

Features and benefits

- Chemically & Biologically inert
- High Temperature Resistant (500°C)
- High flow rate & high dirt holding capacity

Applications

Engineered for air sampling in aggressive environments, the reinforced membrane delivers superior mechanical stability and chemical resistance, making it well-suited for applications involving corrosive aerosols or other demanding testing conditions.

Borosil Filter type Cat no FPBGFEP2000


High-purity micro glass fiber filters are made entirely from borosilicate glass and are specifically designed for use in high-volume PM-10 air sampling systems. These filters effectively capture airborne particles and aerosols, allowing for precise chemical analysis of trace pollutants with minimal interference

BOROSIL PTFE Filter Type-Cat FPBPTPM25

This advanced thin PTFE membrane, known for its high purity and strong chemical resistance, is mounted on a sequentially numbered polypropylene ring. It is perfectly suited for PM2.5 ambient air monitoring and sampling, even in demanding environmental conditions.



The benchmark in Filter Papers

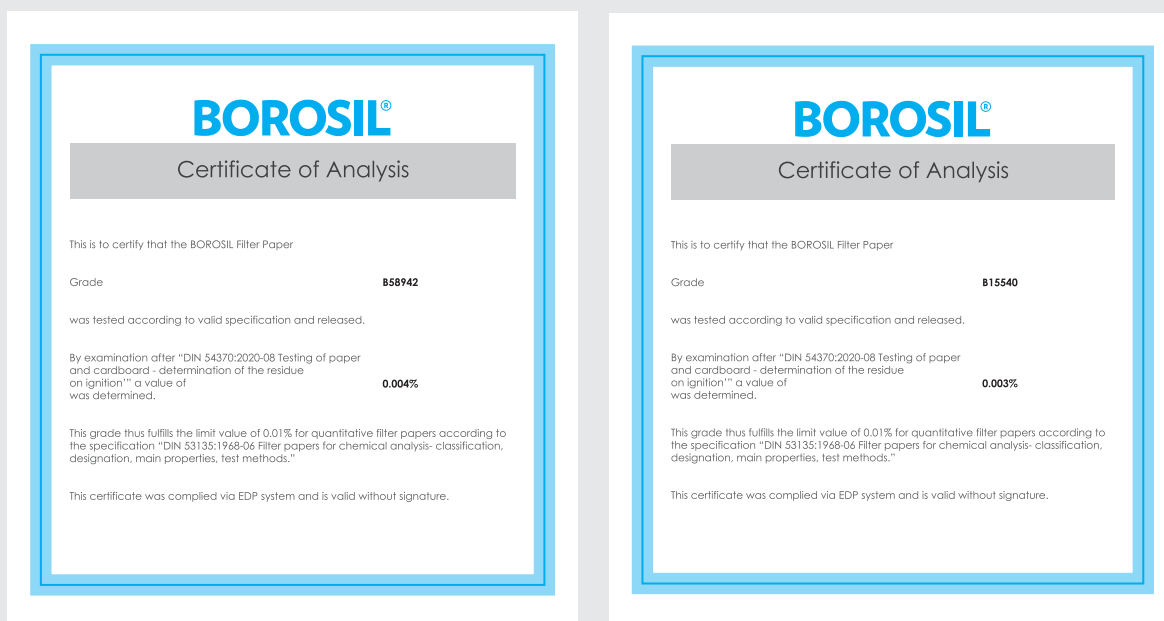

The average ash content of the Quantitative papers is **0.004%** and that for Hardened Ashless is **0.003%**




Applications

- Water pollution analysis: Determination of suspended particles
- Biochemical issues like DNA, RNA, proteins and polysaccharides
- Determination of suspended particles (SPM and TSP)
- Cleaning and buffering solutions and reagents for spectrophotometric measurements
- Measurement of nitrogen (gravimetrically) in emission analysis
- Measurement of metals, inorganic lead in emission control analysis

Ordering Specification

Grade	Size	Pack Size	Product Code
BGF50A	47mm diameter	100	FPBGF50A0047
BGF50A	203*254mm sheet (8*10 inches)	100	FPBGF50A203254
BGF52C	47mm diameter	100	FPBGF52C0047
BGF52C	125mm diameter	100	FPBGF52C0125

Industrial Applications

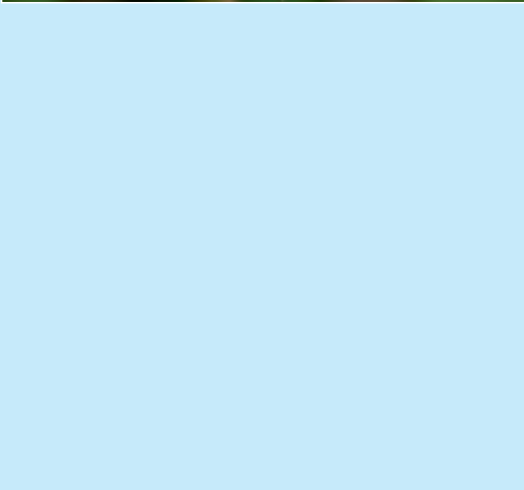

Agriculture



The optimum quality of the soil, fertilizers and animal feed is crucial for the appropriate growth and nutrition of plants and animals. The determination and quantification of nutrients and trace elements hence becomes very critical. The average ash content of BOROSIL quantitative papers therefore make way for accurate, reproducible, error-free results.

Target applications include :

- Analysis of nutrients, mineral nutrients, contaminants and microbiological purity
- Ideal for the detection of trace elements like Mg, Mn, Zn, Co, Cu, Mo, and B

Used in the analysis of :

- Soil and Fertilizers
- Animal feed

	Parameter / Process	Type of Filter	Filter Grade
Soil and fertilizers	Measurement of nitrogen	Measurement of nitrogen	B58940
	Measurement of trace elements	Filter paper for quantitative analysis, ash-free	B58941, B58943, B58942, B58944, B58940
	Free amino acids and total amino acids	Filter paper for quantitative analysis, ash-free	B58943, B58940
	Measurement of soluble sulphates water extraction	Filter paper for quantitative analysis, ash-free	B58942
	Determination of K and P	Filter paper for quantitative analysis, ash-free	B58941, B58943, B58942, B58944, B58940
	Measurement of solids in suspension	Glass microfibre filters	BGF52C
Animal feed	Gravimetric analysis	Filter paper for quantitative analysis	B58941, B58943, B58942, B58944, B58940
	Measurement of Calcium	Filter paper for quantitative analysis	B58943
	Separation of solids from suspension ^s	Glass microfibre filters	BGF52C



Environmental Analysis



A high quality and pure filter paper is required to simplify and support contaminants-free sampling of air, water etc. BOROSIL filter papers can be confidently used in areas that are subject to strict official requirements like the DIN, EPA, ASTM, etc.

Target applications include:

- Ambient air monitoring
- Determination of suspended particles and total suspended particles in air.
- Monitoring the presence of pollutants in the air at different measuring points.
- Monitoring of anthropogenic atmospheric emissions (oil refineries, power stations, burning of liquid and solid fuels, cement factories, mining industries, incinerators, iron foundries, grinders, asphalt plants, glass makers, ceramic factories) and at stationary sources.
- Measurement of dust release in workplace and production processes, exhaust fumes from private houses, and newly developed engines (for cars and other vehicles).
- Gravimetric analysis of organic and inorganic contaminants in water and waste water.
- Monitoring microbiological quality of drinking water.
- Determination of total dry residue in water.
- Determination of dissolved organic carbon (DOC) and total organic carbon (TOC) in water.
- Analysis of waste products in the disposal of industrial waste and laboratory waste

The papers are used to analyze :

- Air pollution
- Emission control
- Water
- Waste products

	Parameter / Process	Apparatus	Technique	Type of filter	Filter Grade
Air Pollution	Sampling of total suspended particulate matter TSP ($\varnothing > 30\mu\text{m}$) 1)	High volume capturer		Glass microfibre filters, in line with US EPA	BGF50A
		Low volume capturer			BGF50A
		Cascade impactor			BGF50A
Emission Control	Measurement of nitrogen (gravimetry)	Isokinetic probe with rear filter-holder (up to 500°C)	Filtration, weighing	Glass microfibre filters	BGF50A
	Measurement of inorganic lead	Isokinetic probe with rear filter-holder (up to 500°C)	Atom absorption spectroscopy	Glass microfibre filters	BGF50A
	Measurement of metals	Isokinetic probe with rear filter-holder (up to 500°C)		Glass microfibre filters	BGF50A
	Measurement of deposition of radioactive aerosols	Filtering instrument	Filtration, Scintillation	Glass microfibre filters, retention capability < 1 μm	BGF6
	Monitoring the combustion air	Monitoring the combustion air	Filtration, weighing	Glass microfibre filters	BGF8, BGF9
	Monitoring particles in air and gases	Automatic air filter units, air analysers with filter rolls	Filtration, weighing	Glass microfibre filters with high mechanical strength	BGF10
Water	Sample preparation		Clarification	Qualitative filter paper	B59501
	Total dry residue, ash residue		Filtration, weighing	Glass microfibre	BGF6
				Quantitative filter paper	B58941
	Measurement of solids in suspensions after drying at 105°C		Filtration, weighing	Glass microfibre	BGF52C, BGF6
	Measurement of the total remainder after drying at 180°C			Glass microfibre	BGF52C, BGF6
	Measurement of solids and volatiles after incineration at 550°C			Glass microfibre	GF50A
	Colouration, Radioactivity, Measurement of metals		Filtration	Glass microfibre	BGF6, BGF50A
	Measurement of total and dissolved organic carbon		Filtration, oxidation	Glass microfibre	BGF6, BGF52C
	Measurement of metals				B58941, B58942
Measurement of non metallic inorganic compounds		Filtration	Quantitative filter paper	B58941, B58942, B58940	
Waste Products	Characterisation of toxic substances		Pressure filtration	Glass microfibre filters	BGF52C

Chemicals



Quality is the utmost critical parameter for any chemical reagent, chemicals or pharmaceutical substance. The quality of these chemicals define the quality of the end product and hence the overall success for the company. Superior quality products therefore need accurate testing methods with materials that produce precise, error-free and reproducible results.

The quantitative filter papers of BOROSIL are the purest paper in the filter market. The average ash content is between 0.005%, and 0.003%.

Target applications include :

- Filtration before quantitative analysis
- Sample preparation before HPLC
- Microbiological investigations
- Gravimetric measurements of elements and contaminants.
- Determination of grind level in cement.

Used in the analysis:

- Chemicals
- Detergents
- Oil refinery
- Cement
- Metal ores and minerals

Beverages



To ensure beverages are optimum for consumption in this highly competitive and health conscious world, beverages undergo several QC steps for checking and maintaining appropriate quality of the product.


Target applications include:

- Analysis of ingredients, contaminants and microbiological purity
- Particle separation and clarification before optical measurements
- Sample preparation before sensitive analyses such as HPLC Process filtration
- Measurement of nitrogen compounds, proteins and trace elements in beer, malt and beer based beverages.
- For removing CO₂ and turbidities during beer production.

Used in the analysis of:

- Juice
- Wine
- Beer, wine, and beer based beverages

	Parameter / Process	Technique	Type of filter	Grade
Quality Control	Gravimetry	Filtration (funnel, Büchner)	Filter paper for quantitative analysis	B58941, B58943, B58942, B58940, B58944
			Hardened filter paper for quantitative analysis	B15540, B15542
	Clarification of samples	Pre-filters for membranes	Glass microfibre filters	BGF9
Cleaning Materials	Gravimetry	Filtration (funnel, Büchner)	Filter paper for quantitative analysis	B58941, B58943, B58942, B58940, B58944
	Determination of tenside content	Filtration (Funnel/Büchner)	Glass microfibre filters	BGF50A
	Separation of solids in suspensions	Filtration (Funnel/Büchner)	Glass microfibre filters	BGF52C
Oil Refinery	Gravimetry	Filtration (funnel/Büchner)	Filter paper for quantitative analysis	B58941, B58943, B58942, B58940, B58944
			Hardened filter papers for quantitative analysis	B15540, B15542
Cement Analysis	Blaine test (grinding fineness of cement)		Filter paper for quantitative analysis	B58941, B58943

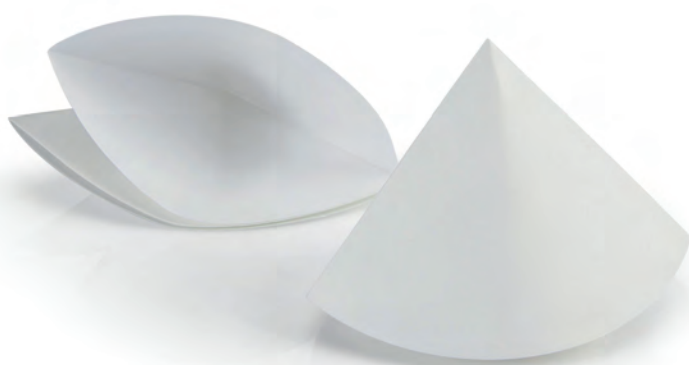

	Parameter / Process	Type of filter	Grade
Juice	Particle separation	Filter paper for qualitative analysis, low ash	B60404, B59702, B59501, B59303
	Preparing fruit juice samples for photometric measurements (e.g. phosphate)	Quantitative filter paper	B58941
	Spectrophotometry clarification of samples	Glass microfibres	BGF6, BGF5
Wine	Particle separation	Filter paper for qualitative analysis	B60404, B59702, B59501, B59303
	Gravimetric analysis	Filter paper for quantitative analysis	B58942
	Sample preparation Pre-filtration	Glass microfibre filters	BGF51B
Beer, malt and beer - based beverages	Removal of CO ₂ and turbidities from beer, wine and juices	Filter paper for qualitative analysis	B59702
	Determination of solids in wort (Labor Veritas method)		BGF52C
	Filtration of lees		B59702, BGF52C
	Determination of the coagulateable proteins		B59702
	Determination of the grade of fermentation		B59702
	Sample preparation		B59501
	Determination of solids and turbidity (Feld method)	Filter paper for quantitative analysis	B58941
	Determination of nitrogen compounds by phosphor molybdenum precipitation		B58943
	Determination of carbohydrates by hydrolysis		B58943
	Analysis of ash content in foodstuffs		B58941
	Determination of proteins in wort and beer via MgSO ₄ precipitation		B58941, B58943
	Drinking water: Determination of chemical elements, radioactive trace elements		B58942





Knowledge about the components and ingredients in various foodstuffs is extremely essential for consumers.

BOROSIL high purity and low ash filter papers are perfectly suitable and compatible with the food analysis tests, and allow detections of sensitive contaminants and ingredients, hence guaranteeing accurate results

Target applications include:

- Analysis of ingredients, contaminants and microbiological purity
- Analysis of oil and components
- Filtration of various extracts
- Determination of fats, proteins, added sugars in milk and milk products

Used in the analysis of

- Edible oil
- Sugar
- Milk and milk products
- Meat and meat products

	Parameter / Process	Type of filter	Grade
Edible oil and fat	Determination of the unsaponifiable fraction in fat	Filter paper for qualitative analysis	B59702, B59501
	Analysis of oil/fats using fat extracting equipment	Filter paper for quantitative analysis	B58940
Sugar	Gravimetric analysis	Filter paper, quantitative analysis	B58941, B58943
	<i>Improvement in filtration-Clarification of samples-Pre filters for membranes</i>	Glass microfiber filters	BGF9
Milk and milk products	Qualitative Analysis	Filter paper for qualitative analysis	B60404, B59501, B59702
	Gravimetric analysis	Filter paper for quantitative analysis	B58941, B58943, B58942
	Measurement of solids in suspensions	Glass microfibre filters	BGF52C
Meat and meat products	Gravimetric analysis	Filter paper for quantitative analysis	B58941, B58943, B58942, B58940



Pharmaceuticals



Pharmaceutical and Healthcare products require stringent QC and papers that do not interfere with the analysis of substances. The low ash content of BOROSIL quantitative papers ensure error-free results along with consistent filtering performance.

Target applications include :




- Monitoring purity, contamination and inspecting microbiological purity
- Clarification before analysis
- Gravimetric measurements
- Sample preparations before HPLC

Used in :

- Production and analysis of APIs
- Quality control of final drug products, healthcare products
- Analysis of various excipients used during production of healthcare and cosmeceuticals

	Parameter / Process	Technique	Type of filter	Grade
Production and Quality Control	Gravimetry	Filtration (funnel, Büchner)	Filter paper for quantitative analysis	B58941, B58943, B58942, B58940, B58944
	Clarification of samples	Pre-filters for membranes	Pre-filters for membranes	BGF9
Diagnostics	Separation and isolation of DNA, RNA	Filtration (even by centrifuge)	Glass microfibre without binder	BGF50A, BGF51B, BGF52C
	Tests for diseases and allergies etc	Sample device for detection reactions with enzymes, antibodies (impregnation)	Pure, absorptive filter papers, almost no contaminants	B58941, B58943, B58942, B58940, B58944



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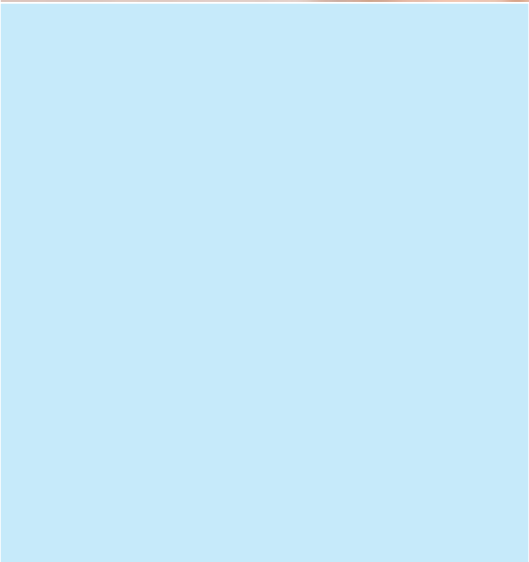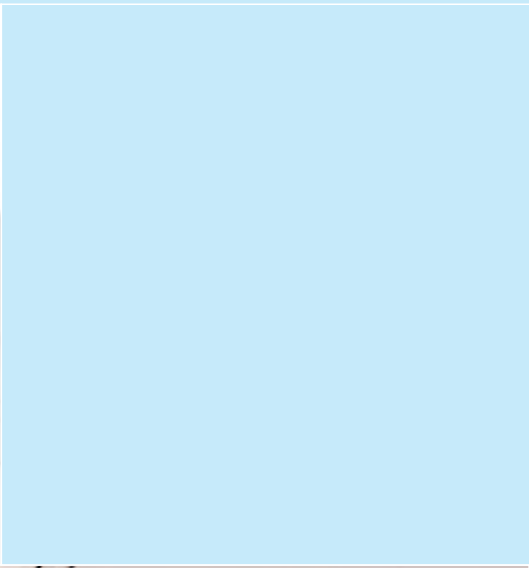



CIN : U74999MH1991PLC061851

Visit
Website



Follow us on
LinkedIn



Registered & Corporate Office :

1101, Crescenzo, G-Block, Opp. MCA Club, Bandra Kurla Complex,
Bandra (E), Mumbai - 400 051, India.

T +91 22 6740 6300 ● F +91 22 6740 6514

E bsl@borosil.com ● W www.borosilscientific.com

Write to us for export enquiries at bsl.exports@borosil.com

MUMBAI

mumbaisales@borosil.com

+91 22 6740 6400

+91 22 6740 6444

DELHI

delhi@borosil.com

+91 11 4150 5893

CHENNAI

chennai@borosil.com

+91 44 2822 6012 / 13

+91 44 2822 6014

KOLKATA

calcutta@borosil.com

+91 33 2229 9166

All rights reserved. Borosil Scientific Limited. 2025.

Any unauthorized use, copying, or distribution of this brochure is strictly prohibited and constitutes a legal violation. Borosil Scientific Limited reserves all rights to take legal action against individuals or entities who violate these legal rights.